



WF300 Quick Installation instructions

1. According to the size of the microscope eyepiece diameter, select the appropriate adapter(seven adapters, diameter from 20 ~ 46mm); for eyepiece diameter in 42-46mm, no adapter is needed.


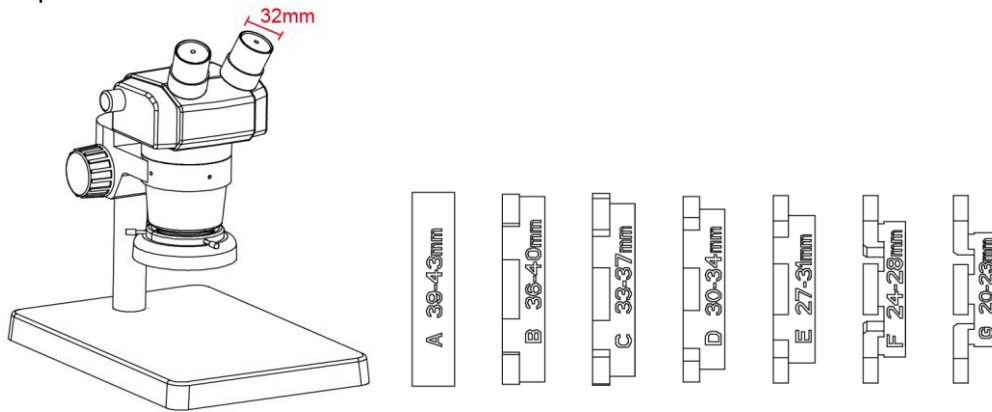


2. For example: eyepiece diameter , choose adapter # D (30-34mm) , put into the chuck, remove the eyepiece, flat on the inner jacket,tighten the swivel.

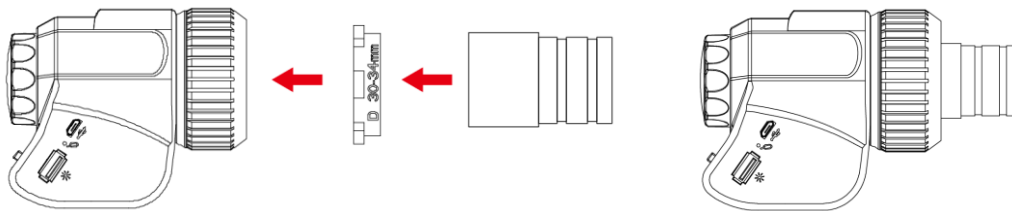

3. Insert into microscope eyepiece tube, start work modes(both USB and WiFi work modes available).

4. Adjust the focal length of the exit pupil until the tangent circle white image Consistent with the visual effect. Adjust the microscope working distance and zoom knob to the proper position. Installation complete.

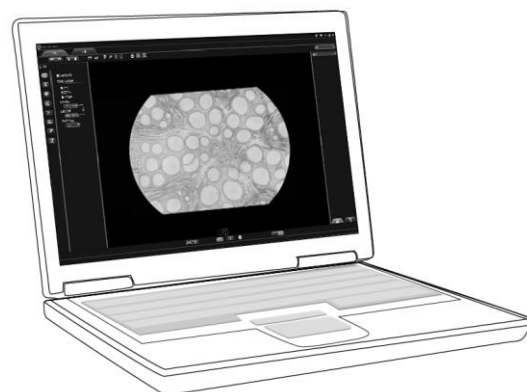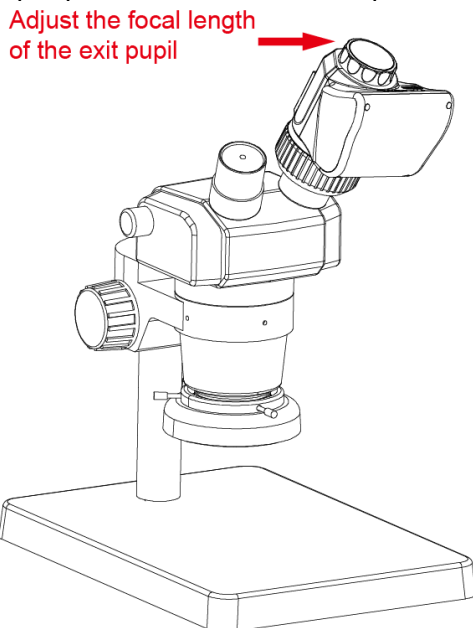

5895 Rue Ferrari San Jose, CA 95138
1 (800) 832-0060 | Fax: (408) 226-0900
www.meijitechno.com